This simplified safety manual describes only the items necessary for using this product safely. You can obtain the SZX-AR1 system instruction manual that describes the details of setup procedures or specifications, etc. from the "Instruction Manual Downloads" website shown below. Before using this product, download the instruction manual from the website and read it carefully.

To read the instruction manual, it is necessary to install software for browsing the PDF file on the terminal you use in advance. In addition, if you cannot browse the instruction manual, do not use the product.

Instruction Manual Downloads

<https://www.olympus-ims.com/service-and-support/instruction-manual-downloads/>

Service Center

<https://www.olympus-ims.com/service-and-support/service-centers/>

Our Website

<https://www.olympus-ims.com>

The following symbol is used in this manual.

⚠ WARNING:

Indicates a potentially hazardous situation which, if not avoided, could result in death or serious injury.

⚠ CAUTION:

Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices or potential material damage.

NOTE:

Indicates a potential situation which, if not avoided, may result in failure of this instrument.

⚠ WARNING: Prevention of electric shock

Never disassemble any part of this product.

It could cause electric shock or failure of the product.

Do not touch the product with wet hands.

In particular, if you touch the main switch of the power unit or the power cord with wet hands, electric shock, ignition or failure of the product may be caused.

Do not bend, pull or tie the power cord/cables in a bundle.

Otherwise they could be damaged, causing a fire or an electric shock.

Keep the power cord and cables well away from the lamp housing.

If the power cord and cables contact a hot area of the lamp housing, they could melt and cause electric shock.

⚠ CAUTION: Electric safety

Always use the power cord specified by us.

If the proper AC adapter and the power cord are not used, the electric safety and the EMC (Electromagnetic Compatibility) performance of the product can not be assured. If no power cord is provided, please select the proper power cord by referring to the section "Proper selection of the power cord" at the end of the instruction manual (Hardware).

Always connect the ground terminal.

Connect the ground terminal of the power cord and that of the power outlet.

If the product is not grounded, our intended electric safety and EMC performance of the product can not be assured.

Do not use the product in close proximity to the sources of strong electromagnetic radiation.

Proper operation may be interfered. The electromagnetic environment should be evaluated prior to operation of the product.

Disconnect the power cord in case of emergency.

In case of emergency, disconnect the power cord from the power cord connector on the product or from the power outlet.

Install the product at the location where you can reach the power cord connector or the power outlet at hand to disconnect the power cord quickly.

Do not connect or disconnect the power cord, cables and units while the power is on.

Safety symbols


The following symbols are placed on this product.

Study the meaning of the symbol and always use the product in the safest possible manner.

Symbol	Meaning
	Indicates a non-specific general hazard. Follow the description given after this symbol or in the instruction manual
	Indicates that the seesaw type main switch is ON. (Seesaw type is the type of switch where ON or OFF is selected by pressing it to ON or OFF side)
	Indicates that the seesaw type main switch is OFF.

Caution engraving/label

Safety labels are placed on the portions which particularly require special cautions when using or operating this product. Be sure to follow these instructions.

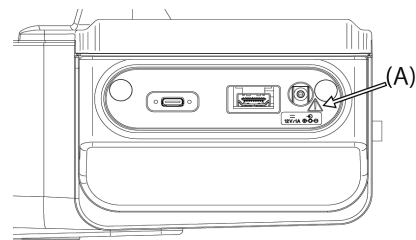Label position	Safety label	Instructions in the instruction manual
(A)		[Caution for electric safety] [Caution for prevention of electric shock]

If a caution engraving or label is dirty or peeled off, contact us for the replacement or other inquiry.

Maintenance and Storage

Cleaning of each part

Lens and filter

Do not leave stains or fingerprints on the lenses or filters. If they get dirty, blow away dust with a commercially available blower and gently wipe the lens or filter with a piece of cleaning paper (or washed-out clean gauze). Only when cleaning fingerprints and oil stains, slightly moisten a piece of cleaning paper with commercially available absolute alcohol and wipe them off with it.

⚠ WARNING:

Since the absolute alcohol is highly flammable, it must be handled carefully. Be sure to keep it away from open flames or potential sources of electrical sparks.
For example, the electrical equipment that is switched ON and OFF may cause the ignition of a fire.
Also, always use the absolute alcohol only in a well-ventilated room.

Specifications

Rating

AC adapter	Input : 100-240V \sim 47-63Hz 0.93A(Max) Output: 12.0V --- 3.33A(Max)
AR Tilting Trinocular Tube	Input : 12V --- 1A(Max)

Operating environment

Temperature	10 to 35 °C
Humidity	20 to 85%
Variation of power supply voltage	Rated voltage: \pm 10 %
Pollution degree	2 (in accordance with IEC60664)
Installation category (excess voltage category)	II (in accordance with IEC60664)
Installation location	Tilt of stand: \pm 3° or less
Conditions for safety standards	<ul style="list-style-type: none">• Indoor use• Altitude: Max. 2000 meters• Temperature: 5 to 40 °C• Relative humidity: 20 to 80% for temperatures up to 31 °C (without condensation) In case of over 31 °C, the relative humidity is decreased linearly through 70% at 34 °C, 60% at 37 °C, and to 50% at 40 °C.

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